





HighFidelity Cy3 PCR Labeling Kit

Preparation of Cy3-labeled DNA probes by PCR

Cat. No.	Amount
APP-101-CY3-S	10 reactions x 20 µl
APP-101-CY3-L	50 reactions x 20 µl

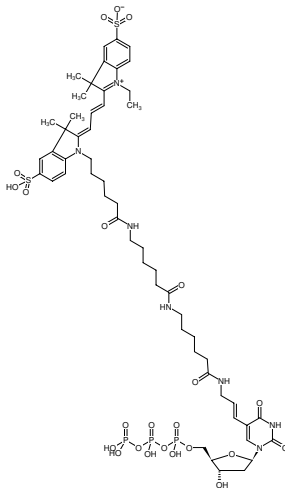

Structural formula of HighFidelity Cy3 PCR Labeling Kit

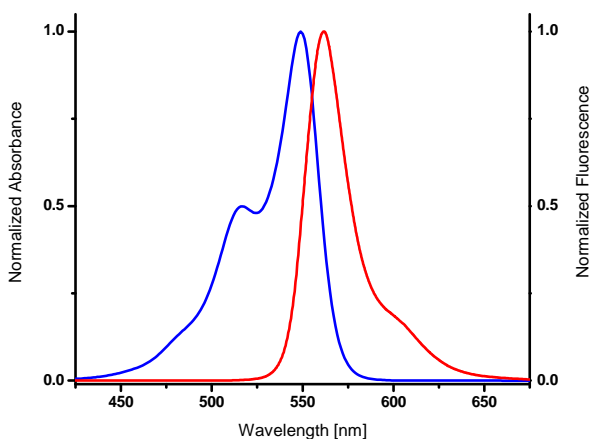

excitation and emission spectrum of Cy3

For general laboratory use.

Shipping: shipped on gel packs

Storage Conditions: store at -20 °C

Additional Storage Conditions: avoid freeze/thaw cycles, store dark

Shelf Life: 12 months

Spectroscopic Properties: λ_{exc} 550 nm, λ_{em} 570 nm, ϵ 150.0 L mmol⁻¹ cm⁻¹ (Tris-HCl pH 7.5)

Description:

HighFidelity Cy3 PCR Labeling Kit is designed to produce randomly Cy3-modified DNA probes by PCR. Such probes are ideally suited for Fluorescence *in situ* hybridization (FISH) and Northern Blot experiments. PCR-based labeling is superior to random-primed labeling with Klenow fragment if template amounts are limited or amplification of a specific DNA fragments is required. Amplification of probes up to 4kbp is feasible.

dUTP-XX-Cy3 is efficiently incorporated into DNA as substitute for its natural counterpart dTTP using an optimized reaction buffer and a High Fidelity Polymerase blend consisting of *Taq* polymerase and a proofreading enzyme. 50 % dUTP-XX-Cy3 substitution typically results in an optimal balance between reaction and labeling efficiency. Individual optimization of dUTP-XX-Cy3/dTTP ratio however, can easily be achieved with the single nucleotide format.

The kit contains sufficient reagents for 10 labeling reactions (S-Pack) or 50 labeling reactions (L-Pack) of 20 µl each (50% dUTP-XX-Cy3 substitution, 100 µM dATP/dGTP/dCTP, 50 µM dTTP, 50 µM dUTP-XX-Cy3).

Content:

High Fidelity Polymerase

in storage buffer with 50% glycerol (v/v)

#APP-101-Cy3-S: 1x 40 µl (100 units, 2.5 units/µl)

#APP-101-Cy3-L: 2x 40 µl (2x 100 units, 2.5 units/µl)

High Fidelity Labeling Buffer

1x 500 µl (10x)

dATP - Solution

1x 20 µl (100 mM)

dGTP - Solution

1x 20 µl (100 mM)

dCTP - Solution

1x 20 µl (100 mM)



HighFidelity Cy3 PCR Labeling Kit

Preparation of Cy3-labeled DNA probes by PCR

dTTP - Solution

1x 20 µl (100 mM)

dUTP-XX-Cy3

#APP-101-Cy3-S: 1x 10 µl (1 mM)

#APP-101-Cy3-L: 5x 10 µl (1 mM)

Lambda DNA

1x 20 µl (100 ng/µl)

500 bp forward primer

1x 20 µl (10 µM)

500 bp reverse primer

1x 20 µl (10 µM)

PCR-grade water

1x 1.2 ml

To be provided by user

DNA template

Primer

DNA purification tools (optional)

1. Preparation of working solutions

1.1 Preparation of 1 mM dATP/dCTP/dGTP working solution

- Thaw 100 mM dATP, 100 mM dCTP and 100 mM dGTP solutions on ice, vortex and spin-down briefly.
- Prepare a 1:100 dilution with PCR-grade water to achieve a final concentration of 1 mM (e.g. 2 µl 100 mM dATP + 2 µl 100 mM dCTP + 2 µl 100 mM dGTP + 194 µl PCR-grade water).
- 1 mM ATP/CTP/GTP working solution can be stored at -20°C. Prepare aliquots to avoid freeze/thaw cycles.

1.2 Preparation of 1 mM dTTP working solution

- Thaw 100 mM dTTP solution on ice, vortex and spin-down briefly.
- Prepare a 1:100 dilution with PCR-grade water to achieve a final concentration of 1 mM (e.g. 2 µl 100 mM dTTP + 198 µl PCR-grade water).
- 1 mM dTTP working solution can be stored at -20 °C. Prepare aliquots to avoid freeze/thaw cycles.

3. Standard PCR Labeling protocol

The standard protocol is set-up for labeling of a 500 bp DNA fragment. An optimal balance between reaction and labeling efficiency is typically achieved with 50% dUTP-XX-Cy3 substitution

following the standard protocol below however, individual optimization might improve results for individual applications.

- Assemble the PCR on ice in the order stated below (DNase-free reaction tube).
- Vortex and spin-down briefly.
- Perform assay set-up and reaction under low-light conditions.

Component	Volume	Final concentration
PCR-grade water	X µl	
High Fidelity Labeling Buffer (10x)	2 µl	1x
1 mM dATP/dCTP/dGTP working solution (s. 1.1)	2 µl	100 µM
1 mM dTTP working solution (s. 1.2)	1 µl	50 µM
1 mM dUTP-XX-Cy3	1 µl	50 µM
forward primer (10 µM)	X µl	0.1 - 1 µM (e.g. 0.3 µM 500 bp forward primer)
reverse primer (10 µM)	X µl	0.1 - 1 µM (e.g. 0.3 µM 500 bp reverse primer)
template DNA	X µl	1 - 10 ng genomic DNA (e.g. 1 ng Lambda DNA)
High Fidelity Polymerase (2.5 units/µl)	1 µl	2.5 units
Total volume	20 µl	

Recommended cycling conditions

Cycle step	Temperature	Time	Cycles
Initial denaturation	95°C	2 min	1x
Denaturation	95°C	20 sec	30x
Annealing ¹⁾	58°C	30 sec	
Elongation ²⁾	68°C	60 sec	
Final Elongation	68°C	2 min	1x

¹⁾The annealing temperature depends on the melting temperature of primers used.

²⁾The elongation time depends on the length of fragments to be amplified. A time of 2 min/kbp is recommended. Elongation at 72°C



HighFidelity Cy3 PCR Labeling Kit

Preparation of Cy3-labeled DNA probes by PCR

works as well.

For optimal amplification results and high incorporation rates an individual optimization of the recommended PCR assay and cycling conditions may be necessary for each new primer-template pair.

4. Probe purification:

Probe purification is not required for most hybridization experiments. If a downstream application requires purification (e.g. concentration determination by absorbance measurement) we recommend silica-membrane or gel filtration-based purification.

Related Products:

Aminoallyl-dUTP-XX-Cy3, #NU-803-XX-Cy3